

QUICK-TAP – FIXATION SYSTEM FOR PLAIN BACKED PAPER

■ Check the flatness of your working disc and ensure that Quick-Tap is centered properly on the magnetic foil.

■ Remove the release liner from the adhesive layer before applying the plain backed paper.

1 Always position the sample holder so that the samples extend beyond the disc's edge during grinding and fill the sample holder symmetrically.

2 Cleaning: If dirty or decreasing adhesion - use warm water and a few drops of dish soap together with a soft sponge.

3 Rinse with clean water and dry by boost function or air.

■ To avoid dust on the adhesive layer, cover the Quick-Tap surface after use with unused plain backed paper. Do not leave uncovered.

